

PPH33 PODS[®] Human NT-4

Description

The product contains the polyhedrin protein co-crystallized with Human NT-4. Neurotrophin-4 (NT-4), also known as NT-5, is a neurotrophic factor and a member of the NGF family. NT-4 is expressed in the prostate, thymus, placenta, and skeletal muscle. NT-4 promotes the development and survival of peripheral and CNS neurons and the proliferation of keratinocyte.

Length	176 aa
Molecular Weight	38.4 kDa
Source	<i>Spodoptera frugiperda (Sf9) cell culture</i>
Accession Number	P34130

Usage Recommendation

PODS[®] co-crystals provide a depot of proteins which are steadily secreted. It has been estimated that the biological activity of 50 million PODS[®] co-crystals generates the same peak dose as 3.3 µg of standard recombinant protein. However, at 5 days following the start of seeding the PODS[®] co-crystals, there are more than 50% of these peak levels still present in the culture system. Ultimately, the amount of PODS[®] co-crystals that is optimal for a particular experiment should be determined empirically. Based on previous data, we suggest using 50 million PODS[®] co-crystals in place of 3.3 µg of standard growth factor as a starting point. To control for cross-reactivity with cells or as a negative control, we recommend using PODS[®] growth factors alongside [PODS[®] Empty crystals](http://www.cellgs.com/products/podsand8482-empty.html), as the latter do not contain or release cargo protein.

Specifications

Alternative Names	Neurotrophin 4, neurotrophin-4, NT4, GLC10, neurotrophic 5, neurotrophin-5, neurotrophic factor 5, neurotrophic factor 4, NT-4/5, NT-5, NT5, NTF4, NTF5
Endotoxin Level	<0.06 EU/ml as measured by gel clot LAL assay
Formulation	PODS [®] were lyophilized from a volatile solution
AA Sequence	MADVAGTSNR DFRGREQRLF NSEQYNNNS KNSRPSTSLY KKAGFMGVSE TAPASRRGEL AVCDAVSGWV TDRRTAVDLR GREVEVLGEV PAAGGSPLRQ YFFETRCKAD NAEEGGPGAG GGGCRGVDRR HWVSECKAKQ SYVRALTADA QGRVGRWIR IDTACVCTLL SRTGRA

Preparation and Storage

Reconstitution	PODS [®] co-crystals may be reconstituted at 200 million co-crystals/ml in water. 20% glucose has a buoyant density closer to PODS [®] co-crystals and can be useful for aliquoting. PODS [®] co-crystals are highly stable when stored in aqueous solution (pH range 6 - 8).
Stability and Storage	Upon receipt, store at 4°C. PODS [®] co-crystals are stable for at least 1 year when dry and 6 months when resuspended.